

Instructions for installing the QX3 Microscope with Windows XP or Vista

Note: DO NOT use the software from the CD that came with the microscope.

DO NOT plug in the QX3 microscope until Step 4.

Step 1

Install the QX3+ download from Intel:

http://downloadcenter.intel.com/Detail_Desc.aspx?agr=Y&Inst=Yes&ProductID=160&DwnldID=4419&strOSs=All&OSFullName=All%20Operating%20Systems&lang=eng

Step 2

Download and install the QCODINSTL Codecs software from:

http://downloadcenter.intel.com/Detail_Desc.aspx?ProductID=160&DwnldID=2089&lang=eng

Step 3

Download and extract the QX3 Drivers from:

<http://hyperslug.s3.amazonaws.com/IntelQX3driverXP.zip>

Step 4

Now, plug in the microscope.

When the "Found New Hardware" wizard appears, select "Locate and install driver software (Recommended)", a message "Installing device driver software" will appear. If you click on "details", you will notice see "Searching Windows Update", after a while you will be able to select "I don't have a disc. Show me other options" and browse to the folder with the extracted QX3 drivers from Step 3. above. The hardware should install successfully.

Double-click on the QX3+ icon on the desktop and the program should run.